



The automatic knife sharpener Leica SP9000 is a highly efficient instrument for resharpening microtome knives, giving a used knife edge the sharpness of a new or factory reconditioned knife. For resharpening, the knife is clamped in the knife holder at a fixed angle. The transparent plastic cover of the instrument is closed during the honing procedure to ensure safe operation and to avoid that dust and grinding particles are scattered over the work place.

Leica SP9000

Automatic knife sharpener

for resharpening

of microtome knives

Leica

MICROSYSTEMS

Leica SP9000 is capable of resharpening all C- profile steel knives up to 25 cm length



Apply a small amount of abrasive to the hone plate as shown in the photograph. Two different types of abrasives are used for the two honing techniques, coarse and fine honing procedures, that are possible.



Place the knife in the knife holder and secure by tightening the clamp screws. Close the plastic cover, select the required mode of operation – coarse or fine honing – and set the automatic timer.



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
The automatic knife sharpener first hones one side of the cutting facet, then automatically turns the knife and hones the other side of cutting edge. This cycle is repeated continuously until the time set with the automatic timer has run out. At the end of the time cycle, the knife holder will stop in a raised horizontal position and the knife can be removed conveniently.

Technical specification subject to change.

Up-to-date development, production and quality control procedures certified under DIN EN ISO 9001 ensure highest quality and reliability.

For further information please contact your local Leica sales company

Australia:	Gladesville/NSW	Phone +61 2 9897 9700	Fax +61 2 9817 8358
Austria:	Vienna	Phone +43 1 486 80 50	Fax +43 1 486 80 50 30
Canada:	Richmond Hill/Ontario	Phone +1 905 762 2000	Fax +1 905 762 8937
Denmark:	Herlev	Phone +45 4454 0101	Fax +45 4454 0111
France:	Rueil-Malmaison Cedex	Phone +33 1 4732 8585	Fax +33 1 4732 85 86
Germany:	Bensheim	Phone +49 6251 136 0	Fax +49 6251 136 155
Italy:	Milan	Phone +39 0257 486.1	Fax +39 0257 40 3273
Japan:	Tokyo	Phone +81 3 5421 2804	Fax +81 3 5421 2894
Korea:	Seoul	Phone +82 2 514 6543	Fax +82 2 514 6548
Netherlands:	Rijswijk	Phone +31 70 4132130	Fax +31 70 4132139
People's Republic of China:	Hong Kong	Phone +852 2 564 6699	Fax +852 2 564 4163
Portugal:	Lisbon	Phone +351 1 388 9112	Fax +351 1 385 4668
Singapore:	Singapore	Phone +65 6779 7823	Fax +65 6773 0628
Spain:	Barcelona	Phone +34 93 494 9530	Fax +34 93 494 9532
Sweden:	Sollentuna	Phone +46 8 6254 545	Fax +46 8 6254 510
Switzerland:	Glattbrugg	Phone +41 1 809 34 34	Fax +41 1 809 34 44
United Kingdom:	Milton Keynes	Phone +44 1 908 246246	Fax +44 1 908 609992
USA:	Bannockburn/Illinois	Phone +1 847 405 0123	Fax +1 847 405 0164



...and more than 100 national distributors.

 www.histo-solutions.com

Leica
MICROSYSTEMS

Order No. 0703-1-0-100
©Leica Microsystems Nussloch GmbH, Heideberger Str. 17-19, D-69226 Nussloch,
Phone (06224) 143-0, Fax (06224) 143-200

11/2005